

Genotyping with the 96.96 IFC Using Standard TaqMan Assays

For more information, see the SNP Genotyping Analysis User Guide (PN 68000098) and the Juno System User Guide (PN 100-7070).

Choose a Juno/IFC Controller HX Workflow

Prime	Load and thermal-cycle (PCR)	Image
Juno™	Juno one-step loading and PCR	Biomark™ HD/Biomark or EP1™

Prime	Load	Thermal-cycle (PCR)	Image
Juno or HX	Juno or HX	Juno or FC1™ cyclers	Biomark HD/Biomark or EP1

Prime	Load	Thermal-cycle (PCR) and image
Juno or HX	Juno or HX	Biomark HD/Biomark

Prime the 96.96 IFC

IMPORTANT

- Use the 96.96 Dynamic Array™ integrated fluidic circuit (IFC) within 24 hours of opening package.
- Due to different accumulator volumes, only use 96.96 syringes with 150 µL of control line fluid.
- Control line fluid on IFC or in the inlets makes IFC unusable.
- Load the IFC within 60 minutes of priming.


- 1 Inject control line fluid into each accumulator on the IFC.
- 2 Remove and discard protective film from bottom of IFC.
- 3 Place the IFC into the instrument and run the prime script:
 - Juno: **Prime 96.96 GT**
 - HX: **Prime (138x)**

Prepare 10X Assays

In a DNA-free hood, prepare aliquots of 10X assays using volumes in the following table. Scale up appropriately for multiple runs.

Component	Vol. per inlet (µL)	Vol. per inlet with overage (µL)	Vol. for 50 µL stock
SNP Genotyping Assay Mix (80X*) (Thermo Fisher Scientific)	0.5	0.625	6.25
2X Assay Loading Reagent (Fluidigm PN 100-7611) ●	2.0	2.5	25.0
ROX™ (50X) (Thermo Fisher Scientific PN 12223-012)	0.2	0.25	2.5
DNA-free water	1.3	1.625	16.25
Total	4.0	5.0	50.0

*If you are using 40X SNP assay, double the volume and reduce the DNA-free water. For other starting concentrations of the SNP assay mix, call Fluidigm technical support.

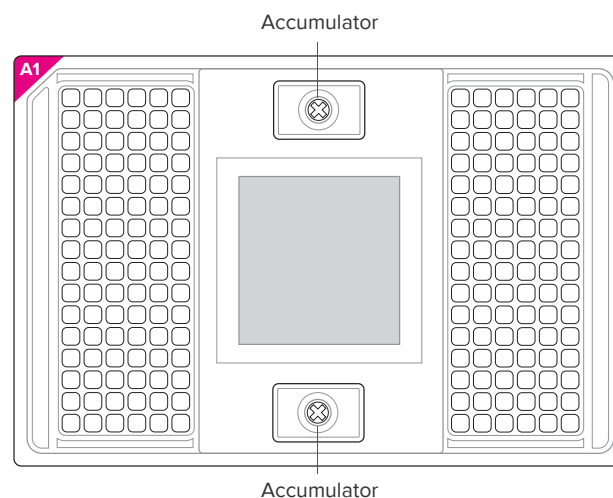

Prepare Sample Pre-Mix and Samples


- 1 Combine components in the following table to make sample pre-mix and final sample mixture.

Component	Vol. per inlet (µL)	Vol. per inlet with overage (µL)	Sample pre-mix for 96.96* (µL)
SAMPLE PRE-MIX			
TaqMan® Universal PCR Master Mix (2X) (Thermo Fisher Scientific PN 4304437)	2.5	3.0	360.0
20X GT Sample Loading Reagent (Fluidigm PN 100-7612) ●	0.25	0.3	36.0
AmpliTaq Gold® DNA Polymerase (Thermo Fisher Scientific PN 4311806)	0.05	0.06	7.2
96.96 GT PCR Dilution Reagent (Fluidigm PN 100-9977)	0.1	0.12	14.4
Genomic DNA (added individually to sample pre-mix)	2.1	2.52	—
Total	5.0	6.0	—

*120 reactions for ease of pipetting

- 2 In a DNA-free hood, combine the four sample pre-mix components in a 1.5 mL sterile tube—enough volume to fill an entire IFC. Aliquot 3.48 µL of the sample pre-mix for each sample.
- 3 Remove the aliquots from the DNA-free hood and add 2.52 µL of genomic DNA to each, making a total volume of 6 µL in each aliquot.

96.96 IFC Pipetting Map

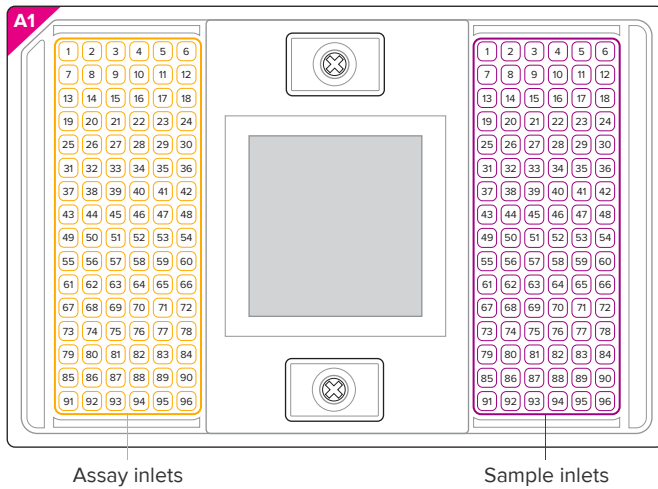

Load the IFC

IMPORTANT

- Vortex thoroughly and centrifuge all assay and sample solutions before pipetting into the IFC inlets. Failure to do so may result in a decrease in data quality.
- While pipetting, do not go past the first stop on the pipette. Doing so may introduce air bubbles into inlets.
- For unused assay inlets, use 2.5 μL assay loading reagent, 0.25 μL ROX and 2.25 μL water per inlet.
- For unused sample inlets, use 3.48 μL of sample mix and 2.52 μL of water per inlet.

- 1 When the prime script has finished, remove the primed IFC from the instrument and pipet 4 μL of each assay and 5 μL of each sample into their respective inlets on the IFC.
- 2 Return the IFC to Juno or HX, then run the script according to operation:

Instrument	Operation	Run script	Continue to
Juno	One-step loading and thermal cycling	One Step 96.96	Collect End-Point Data
Juno	Loading only	Load Mix 96.96 GT	Thermal-Cycle the 96.96 IFC
HX	Loading only	Load Mix (138x)	Thermal-Cycle the 96.96 IFC

IMPORTANT Start IFC run within 1 hour of loading samples.

Thermal-Cycle the 96.96 IFC

Choose the instrument and run the script:

Instrument	Operation	Run script
Juno	One-step loading and PCR	—
Juno	PCR only	Probe GT tab: PCR 96.96
Biomark HD or Biomark	PCR and imaging	Continue to Collect End-Point Data and select GT 96x96 Standard v1.pcl

To run this protocol as an end-point read using the FC1 cyclor or the Fluidigm stand-alone thermal cycler, see the FC1 Cyclor Usage Quick Reference (PN 100-1250) or the Stand-alone Thermal Cyclor Usage Quick Reference (PN 68000111), respectively.

Collect End-Point Data

- 1 Remove any dust particles or debris from the IFC surface.
- 2 Double-click the **Data Collection Software** icon on the desktop to launch the software.
- 3 Click **Start a New Run**.
- 4 Ensure that the status indicators for the lamp (Biomark and EP1 only) and the camera are green.
- 5 Place the IFC into the instrument. Click **Load**.
- 6 Verify IFC barcode and IFC type.
- 7 Choose project settings (if applicable). Click **Next**.
- 8 Provide a name and select a file storage location for a new IFC run, or browse to select a predefined run file. Click **Next**.
- 9 Choose the application, reference, and probes:
 - a Application type: **Genotyping**
 - b Passive reference: **ROX**
 - c Probe types: **FAM-MGB** and **VIC-MGB**
 - d Click **Next**.
- 10 Browse to and choose a thermal protocol:
 - Biomark HD or Biomark for end-point read only (after cycling on Juno or FC1), select **GT End Point v1**
 - Biomark HD or Biomark for thermal cycling and imaging, select **GT 96x96 Standard v1.pcl**
 - Be sure to use a 96.96-specific protocol.
 - EP1, continue to the next step.
- 11 Confirm **Auto Exposure** is selected. Click **Next**.
- 12 Verify the IFC run information. Click **Start Run**.

For technical support visit fluidigm.com/support

North America +1 650 266 6100 | Toll-free: 866 358 4354 in the US | techsupport@fluidigm.com
China (excluding Hong Kong) +86 21 3255 8368 | techsupportchina@fluidigm.com
All other Asian countries +1 650 266 6100 | techsupportasia@fluidigm.com

Europe +33 1 60 92 42 40 | techsupporteurope@fluidigm.com

Japan +81 3 3662 2150 | techsupportjapan@fluidigm.com

Central and South America +1 650 266 6100 | techsupportlatam@fluidigm.com

For Research Use Only. Not for use in diagnostic procedures.

Information in this publication is subject to change without notice. **Safety data sheet information:** fluidigm.com/sds. **Patent and license information:** fluidigm.com/legalnotices. **EU's WEEE directive information:** fluidigm.com/compliance. Fluidigm, the Fluidigm logo, Biomark, Dynamic Array, EP1, FC1, and Juno are trademarks or registered trademarks of Fluidigm Corporation in the United States and/or other countries. All other trademarks are the sole property of their respective owners. © 2016 Fluidigm Corporation. All rights reserved. 06/2016